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Ziegelei 1
D-72336 Balingen
E-Mail: info@kern-sohn.com

Tel: +49-[0]7433- 9933-0
Fax: +49-[0]7433-9933-149
Internet: www.kern-sohn.com

Betriebsanleitung Plattformwaage
Operating instructions Platform balance
Mode d'emploi Balance de plateforme

KERN IOC

Version 1.1
2021-06

Deutsch
English
Français

IOC-BA-def-2111

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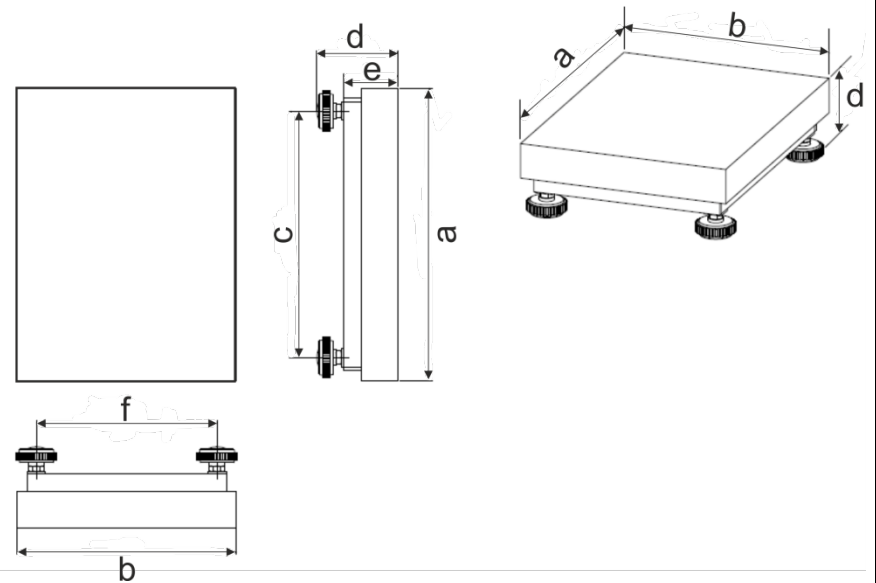
Operating manual Platform

Contents

1	General hints	3
2	Declaration of conformity	3
3	Appliance overview	4
4	Safety instructions	5
5	Assembly of the platform	6
5.1	Levelling	6
6	Cleaning	7
6.1	Packaging / return transport	7
7	Operation	7
8	Cleaning	7
9	Technical data	8
9.1	Dimensions (Measurements in mm)	11

1 General hints

These operating instructions contain all data necessary for placing and commissioning the platform:

The display unit for the platform balance KERN IOC is the model KERN KIB-TM.

Information about

- **Mains connection**
(power is supplied via the connecting cable of the display unit)
- **Initial Commissioning**
- **Connection of peripheral devices**
- **Adjustment and Verification**
(only the complete balance can be verified, i.e. the platform in connection with the display unit KERN KIB-TM).

and the correct operation you will find in the operating instructions included in the scope of delivery of the display unit.

2 Declaration of conformity

The current EC/EU Conformity declaration can be found online in:

www.kern-sohn.com/ce

3 Appliance overview

1. Platform
2. Footscrews
3. Display unit KIB-TM

4 Safety instructions

The product safety has a high ranking at KERN & Sohn.

Non-compliance with the following instructions will result in damage on the platform and / or cause injuries.

- ⇒ Before working on the platform read these instructions.
Keep these instructions for a later use.
- ⇒ Be careful when transporting or lifting heavy devices.
- ⇒ Only qualified personnel may install and care of the platform.
- ⇒ Before cleaning, installation and maintenance, disconnect the platform from the power supply.
- ⇒ The platform must have been stabilized to room temperature, before power supply is switched on.
- ⇒ Do not use the platform in an explosive environment.

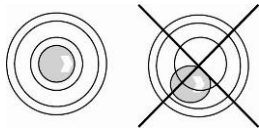
5 Assembly of the platform

On the installation site observe the following:

- Place the platform on a firm, level surface.
The floor at the installation site must be able to carry safely the weight of the maximally loaded platform at the resting points. At the same time it should be sufficiently stable, that no vibrations may occur during weighing work.
- In the installation site possibly no vibrations, e.g. by neighbouring machines should occur.
- Do not use the platform in an explosive environment.
- Avoid extreme heat as well as temperature fluctuation caused by installing next to a radiator or in the direct sunlight.
- Protect the platform against direct draughts e.g. due to open windows and doors.
- Use the platform only in dry environment, protect it against high humidity, vapours and dust.
- Do not expose the device to extreme dampness for longer periods of time. Non-permitted condensation (condensation of air humidity on the appliance) may occur if a cold appliance is taken to a considerably warmer environment. In this case, acclimatize the disconnected appliance for ca. 2 hours at room temperature.
- Avoid jarring during weighing.
- Avoid static charge of goods to be weighed or weighing container.
- Keep away chemicals (such as liquids or gasses), which could attack and damage the balance inside or from outside.
- Keep IP protection of the device.
- Major display deviations (incorrect weighing results) may be experienced should electromagnetic fields (e.g. due to mobile phones or radio equipment), static electricity accumulations or instable power supply occur. Change location or remove source of interference.

5.1 Levelling

Accurate weighing results require a platform with perfect horizontal alignment. During initial installation and after each change of work area it is necessary to level the platform.

- ⇒ As the air bubble is located under the weighing plate, remove it.
- ⇒ Level balance with foot screws until the air bubble of the water balance is in the prescribed circle.

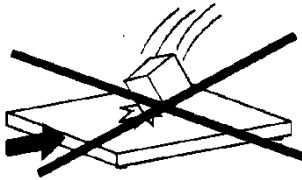
6 Cleaning

6.1 Packaging / return transport

- ⇒ Keep all parts of the original packaging for a possibly required return.
- ⇒ Only use original packaging for returning.
- ⇒ Secure all parts against shifting and damage.

7 Operation

By this way a continuous optimal performance is guaranteed:

- Avoid falling load, shock loads and impacts from the side!
- For weighing operation, all objects must be placed centrally of the weighing bridge and may not hang over the edges.
- Check adjustment at regular intervals.

8 Cleaning

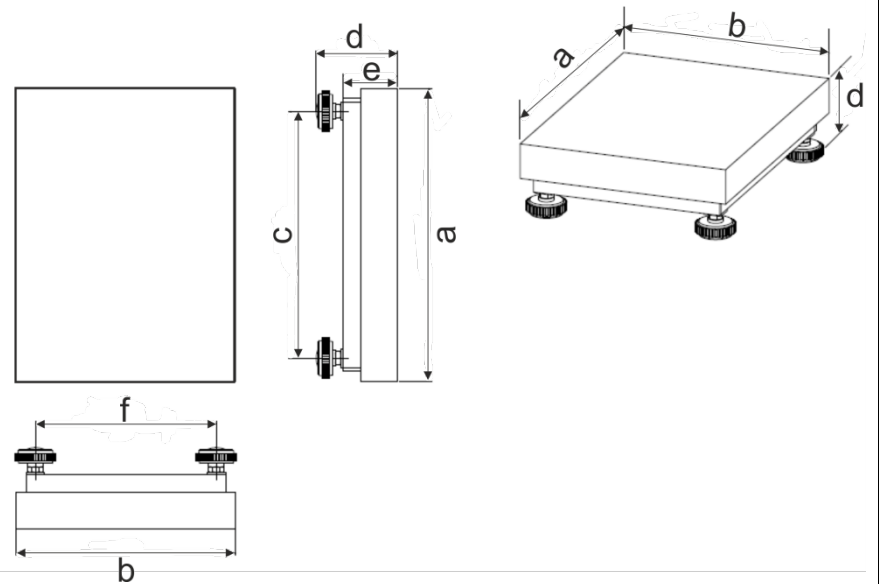
- ⇒ Clean the platform with a soft cloth soaked with a mild cleaning agent.
- ⇒ Take away the weighing plate and remove dirt and foreign matters, which have been accumulated below. Do not use hard or sharp objects for this. Do not open the platform.

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